
IM-55-AV1X FLOW ADJUSTMENT INSTRUCTIONS

Achieve perfectly balanced flow everytime for optimum safety with Method AV1X adjustable eyewash.



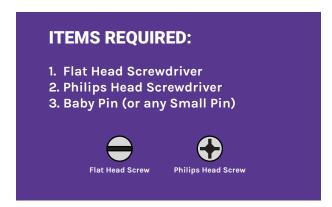

Please follow the below instructions to achieve perfect flow with your AVIX:



Gently remove screw cap with baby pin.



Activate eyewash by pressing down on push handle. Ensure stream of both outlets is balanced (Both stream touching at centre).



Turn gently until desired stream is achieved. Do not force when fully turned. Turn off eyewash and reactivate after 5 seconds. Check result of adjustment.



Use Philips head screw driver to remove regulator lid.



If adjustment needed, adjust regulator by using Flat head screwdriver.



Reinstall regulator lid and screw сар.

3



IM-55-AVIX CLEANING AND SERVICING INSTRUCTIONS

Keep performance at the best level by periodically cleaning your eyewash unit.



ITEMS REQUIRED

- 1. Adjustable Spanner
- 2. Brush with Soft Bristle/ Toothbrush
- 3. Mild Cleaning Soap (Optional)
- 4. Flat Head Screwdriver
- **5.Philips Head Screwdriver**
- 6. Baby Pin (or any Small Pin)



Please follow the below instructions to clean and service your AVIX eyewash:



Gently remove screw cap with baby pin. Use Philips head screw driver to remove regulator lid.



Gently use your thumb to press out honeycomb aerator from the top of eyewash outlet.



Reinstall aerator and eyewash outlet.



Carefully use spanner to detach eyewash outlet.



Rinse aerator under warm water and gently brush until clean. Use soap to remove stubborn grimes.



Finish by performing eyewash stream check. Refer to flow adjustment instructions (2-6).



PERFORMANCE AND ASSURANCE EVERYTIME WITH THE **NEW IM-55-AV1X EYEWASH**

CONVENTIONAL EYEWASH



Occasional unbalanced water stream



Unchecked pressure leads to messy and wet surroundings



Requires supply pressure of 2 bar or more



Parts can be easily tampered or removed

IM-55-AV1X



Consistently achieve perfect stream



Manually regulate flow to ensure steady stream regardless of pressure



Performs well at approximately 1 bar



Uniquely designed to reduce risk of tampering

*IM-55-AV1X eyewash is suitable with previous Method eyewash and shower model.





V1.0-06/21