

FREQUENTLY ASKED QUESTIONS

1) If I have leftover solution can it be kept and reuse in the future?

No, Plum eyewash bottles are for single use only. Once the seal is broken, the risk of contamination makes them unsafe for reuse.

2) Where should I install the station?

- Hazardous work area
- Laboratories
- Manufacturing / Industrial Facilities

Keep in mind install station must be close to high-risk areas (within 10 seconds or 20 meters)*

3) How long should I rinse my eyes?

- For dust or particles: Rinse until discomfort subsides
- For chemical exposure: Rinse continuously for at least 15 minutes and seek medical attention immediately.

4) What is the shelf life of Plum Eyewash?

Shelf life is typically up to 18 months when stored unopened at room temperature. Always check the expiry date printed on the bottle.

5) What is the difference between saline and pH Neutral?

- **Saline (0.9%)** - is ideal for rinsing out dust, dirt, and non-chemical particles.
- **pH Neutral (4.9%)** - is specially formulated to quickly neutralize chemical splashes from acids and alkalis and should be used immediately after chemical exposure.

6) Does Plum Eyewash replace a plumbed emergency eyewash station?


No. Portable Plum Eyewash bottles are intended for immediate first response. In environments with high chemical risk, they should complement, not replace, plumbed eyewash stations.

7) Is Plum Eyewash compliant with safety standards?

Plum products are manufactured in accordance with relevant European safety standards and are suitable for workplace first aid compliance.

8) How to clean Plum Eyewash in case of spill?

In the event of spill, clean it with a damp cloth or mop. No special treatment needed.

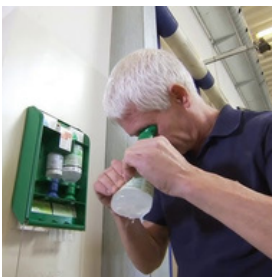

HOW TO USE



1



2



3



4



5

1) Action : Twist the cap to break the seal for immediate use.

2) Positioning : Place the ergonomic eyecup against the eye, covering it completely to hold it open.

3) Application : Squeeze the bottle gently for a soft, continuous flow.

4) Duration : Continue rinsing until the bottle is empty or medical help arrives.

5) Disposal : Plum Eyewash are strictly one time use only. Throw bottle at nearby wastebin.

POPULAR MODELS



EP 3000

Contents:

1x 500ml PLUM Eyewash (EP 4604)
1x Wall Mounted Eyewash Station



EP 4000 + EP 4604

Contents:

1x Wall Mounted Station
2x 500ml PLUM Eyewash (EP 4604)



EP4650

Contents:


1x Dust-proof Wall Station
2x 500ml PLUM Eyewash (EP 4604)



EP 4922

Contents:

1x Portable iBox
2x 500ml PLUM Eyewash (EP 4604)

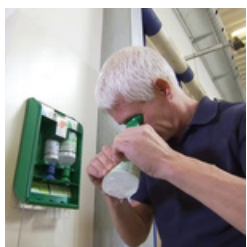

*Contact our salesperson for more information**